

HOW TO SET THE CAMERA ORDER IN VICON NEXUS

Reorder Cameras

1. Turn on the Giganet(s) or Vicon Lock (whatever is applicable to your system).
1. Open Nexus and wait until all devices are booted up properly.
2. Go to the system tab in the Resources window.
3. Right click on 'Vicon Cameras' and select 'Reorder'...

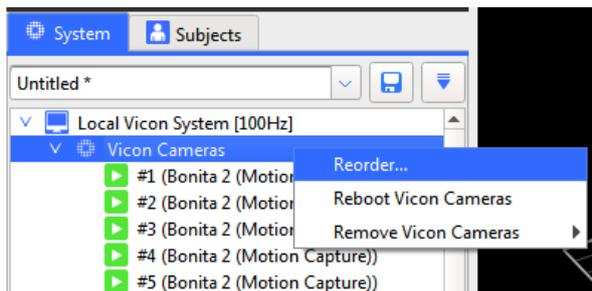


Figure 1: Open the Reorder Devices dialog box

4. The Reorder devices dialog box opens. The cameras are displayed with their current sequence numbers.

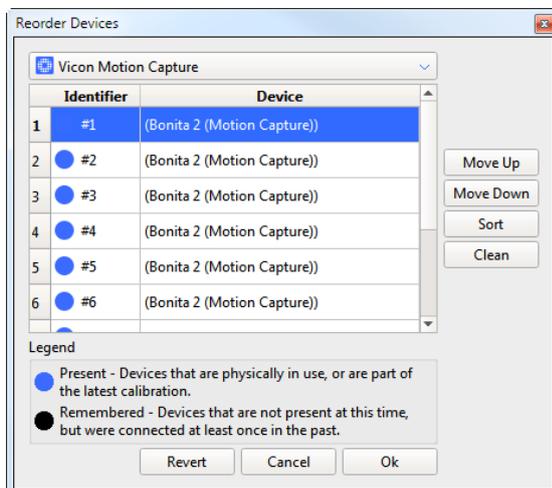


Figure 2: The Reorder Devices dialog box

5. Click on the device, which position you want to change, and move the camera up or down to its desired position. The number of the camera will change immediately.
6. Repeat step 6 for each camera which position you want to change.
7. Click OK to save the changes and close the Reorder Devices dialog box.

Tip: In case the numbering of the cameras does not start with 1, click Sort first.

Tip: First start with the the camera you want to move to position 1. Then find the camera for position 2, and so on.

Determine current camera position

The way to determine the current position is different for Vantage cameras and Bonita cameras.

Determine position of Bonita cameras

1. Bonita cameras do not have an indicator light. To determine the position of a camera set the View pane to 3D perspective, open the Reorder devices dialog box and select a camera in the Reorder devices dialog box.
2. The corresponding camera will be highlighted both in the 3D perspective window as well as in the system Resources pane.

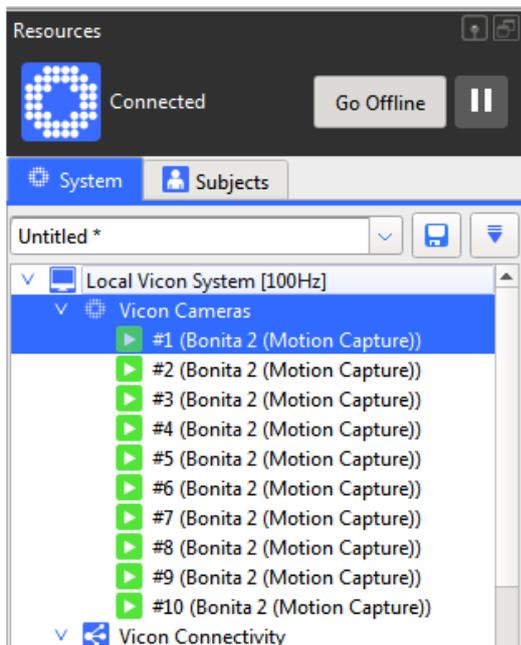


Figure 3: Select a camera in the System Resources pane

Determine position of Vantage cameras

1. In the system resources pane select a camera
2. The indicator light on the camera itself of the corresponding camera will emit a bright, blue light.